




Do-Nature
STV-120M


POCKET MICROSCOPE 60x-120x ZOOM

I) Product overview

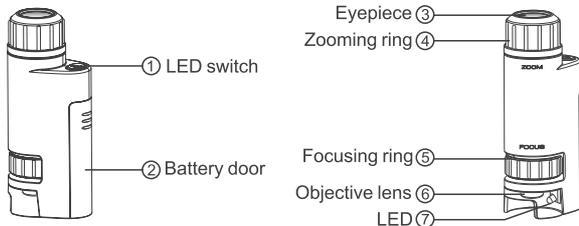

II) Installing the battery


1. Slide the battery door ② as shown below.

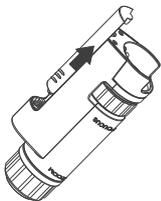

2. Install a AAA battery in the correct polarity.




3. Attach and tighten the battery door ② as shown below.

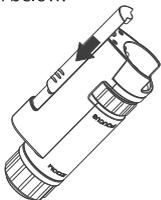

III) Using the microscope

i. Place the viewing object directly below the objective lens ⑥.

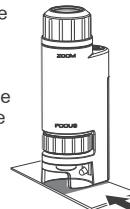

ii. Press the LED switch ① to turn on the LED.



iii. Adjust the zooming ring ④ to the desired level of magnification.



iv. Rotate the focusing ring ⑤ until the image is sharp and clear.



v. Turn off the LED at the end of the observation by pressing the LED switch ①.



IV) Product specifications

Magnification: 60x - 120x

Illumination: LED

Exit pupil diameter: 10mm - 14mm

Size: 37mm x 26mm x 88mm

Optical zoom: 2x

Weight: Approx. 30 grams
(excluding batteries)

Power: 1 x AAA alkaline battery
(not included)

V) Warning

- To avoid possible permanent eye damage, do not look at the sun or bright objects with the microscope.
- To avoid possible permanent eye damage, do not place the lighted microscope directly upon human eye.
- Do not charge non-rechargeable battery; for rechargeable battery, only charge rechargeable battery under adult supervision and remove rechargeable battery from the product before being charged. Remove exhausted battery from the product. Do not short circuit the supply terminals.
- Remove the battery before storing the microscope for an extended period.
- The device may get warm after extended period of use. This is normal and should not be considered as a defect.
- Please retain the packaging for future reference as it contains important information.